


Accurate and Comprehensive Large Molecule Analysis for Quadrupoles

Cerno introduces an entirely new approach for analyzing any large molecule including peptides, oligos, and proteins. SAMMI uses a direct analytical model based on first principles, which out performs the competition by 10-50X. It will analyze each and every ion of a multiply charged molecule to provide not only the neutral mass, but also the abundance distribution across all the charge states.

With SAMMI, you visually observe the results in a simple, intuitive, and confident fashion. Since it is based on a robust and direct analytical model, there are no parameters to tweak, SAMMI is both simple and reliable to use. SAMMI also easily accounts for modifications and adducts and not only determines the accurate neutral mass of the unknown molecule, but also can accurately

confirm the sequence/formula of the molecule as well as its adducts, modifications, impurities, and isotope labels.

SAMMI relies on TrueCal™, Cerno's patented MS calibration method which dramatically improves Spectral Accuracy to allow for the most accurate modeling of the mass spectral profiles for any given ion, regardless of the charge state. It dramatically improves the low resolution MS results to approach the performance from a more expensive high resolution instrument. This makes it perfect not only for research, but for high throughput screening or quality control applications using robust, reliable, and more cost-effective quadrupole instruments. SAMMI works with all large molecules from small peptides and large proteins, oligonucleotides such as RNA of various lengths, and other polymers.

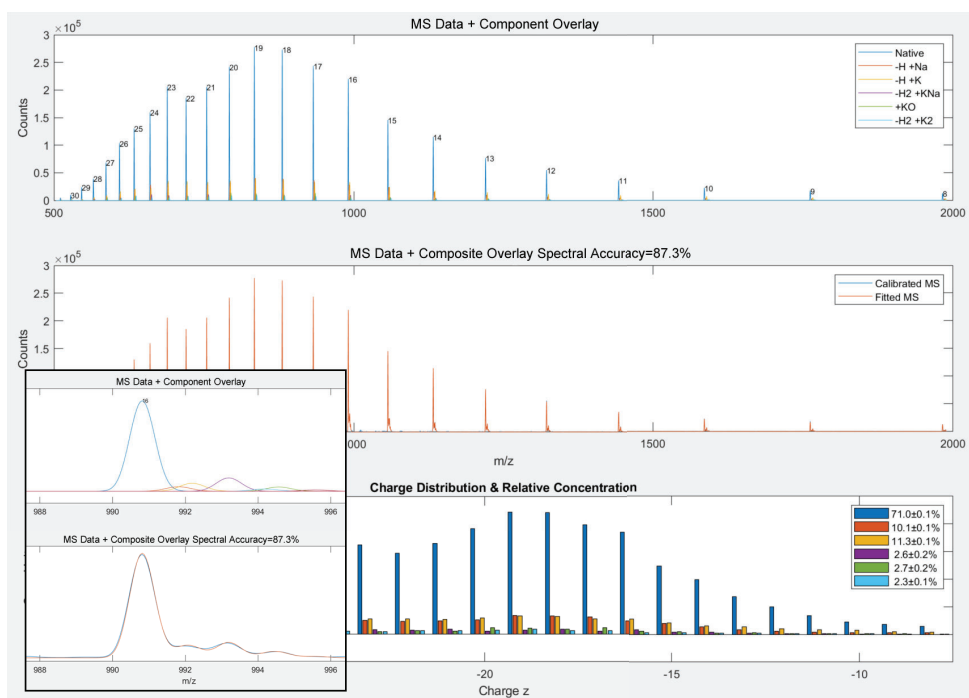

50-nt RNA analyzed against the proposed sequence along with five other components (adducts and modifications, top). The excellent spectral accuracy of almost 90% (middle) confirms the compound, adducts and modifications along with a detailed quantitative relative abundance distribution across all charge states (bottom). Insert shows details for modeled peak at charge state -16.

SAMMI Key Features

- Uses Cerno's proven TrueCal™ approach to directly model large molecules if any charge state
- 10-50X improvement in mass accuracy on single or triple quadrupole instruments
- Accurate mass analysis of unknown molecules as well as their proposed adducts, modifications, and impurities along with all their relative concentrations
- Confirm proposed sequences/formula including multiple modifications and/or isotope labels
- No limit to molecule size, large or small
- Provides relative quantitation for all species across all charge states
- Logical, intuitive, and scientifically defensible interpretation of results, i.e., no more "black box"
- "Parameter Free" making it very easy-to-use
- Available libraries for direct integration with high throughput workflows using a Python API
- Vendor agnostic
- Patented

The Importance of Spectral Accuracy for Large Molecule Analysis

Why is Spectral Accuracy so important to the analysis of large molecules? For the analysis of a singly charged small molecule, a single point measurement of mass accuracy can only confirm a single mass location (either monoisotopic or average mass) among a typically wide mass distribution of the molecules' various isotopes. Spectral Accuracy takes into consideration the entire mass spectral profile including all relevant isotopes and has proven to be a much superior confirmation for the whole molecule. Perhaps more importantly, it can also give a direct indication of any spectral interferences, e.g., the presence of impurities or overlapping ion fragments (e.g. M⁺ and (M-H)⁺ or labeled compounds). With spectra

that are well calibrated with high spectral accuracy, we can explain and quantitatively account for complex multiply charged spectra along with associated adducts, modifications (e.g. deamidation or deamination), and impurities across all charge states. This is accomplished by simply fitting the respective theoretically calculated mass spectra (true MS) to the calibrated mass spectrum through multiple linear least squares regression. Without the use of Spectral Accuracy and by relying solely on mass accuracy, deamidation / deamination or other modifications can easily escape detection, let alone attain accurate and unbiased quantitation, with either low or high resolution instrumentation.

About SAMMI

SAMMI is a companion product to MassWorks, Cerno's award winning calibration and analysis software. Using our new MassLab™ technology, one can add powerful capabilities to MassWorks through custom applications written in Python or MATLAB (MassLab Apps). Currently, SAMMI is a MassLab App available for free with the purchase of or upgrade to the most current version of MassWorks.

cerno
BIOSCIENCE

For More Information

Visit our website at www.cernobioscience.com or email to info@cernobioscience.com
4220 S Maryland Pkwy, Suite 200 · Las Vegas, NV 89119-7523 · USA Tel 1-203-312-1150

